

## (continued) Appendix A. Operator PMCS

4110-01-287-7111  
Refrigerator, Solid State, Biological, Model DLA-50T

[B-Before Operation, D-During Operation, A-After Operation, Q-Quarterly, and S-Semiannually]

ITEM NO	INTERVAL	ITEM TO BE INSPECTED AND PROCEDURE	IS NOT MISSION CAPABLE IF:
1		<b>Refrigerator, Solid State, Biological</b>	
	B, A	a. Conduct an inventory to ensure that the components and accessories listed on the Parts and Accessories List are on hand.	
	B	b. Inspect the unit for dust, dirt, or damage. Inspect power cord for cracks or tears.	The power cord is cracked, frayed, or wires are not covered by the cord insulation.
	B	c. Verify that the door seal is serviceable.	The door does not seal.
	B	d. Verify that the door cover closes and latches properly.	The door cover does not close.
	B	e. Assemble the refrigerator power interconnections as directed in the manufacturer's literature. Ensure that the power source is correct.	
	B, S	f. Install the "Alarm Battery Pack" and test the "Alarm/Battery System" as directed in the manufacturer's literature.	The alarm red light does not blink or the sound signal will not beep.
	B	g. Start the operation of the refrigerator as directed in the manufacturer's literature.	The refrigerator cannot be started.
	B	h. Clean the refrigerator as directed in the manufacturer's literature.	
	A	i. Store the refrigerator as directed in the manufacturer's literature.	