

4110-01-287-7111
Refrigerator, Solid State, Biological, Model DLA-50T

[M-Monthly, Q-Quarterly, S-Semiannually, and A-Annually]

ITEM NO	INTERVAL	ITEM TO BE INSPECTED AND PROCEDURE	IS NOT MISSION CAPABLE IF:
1	S	Refrigerator, Solid State, Biological Conduct an inventory to ensure that the items listed on the Equipment Parts and Accessories List are on hand.	Missing components preclude operation of the refrigerator.
2	S	Maintaining Refrigerator a. Verify that the refrigerator has been maintained according to the Operator Preventive Maintenance Checks and Services. b. Verify that the refrigerator functions on AC current. c. Verify that the refrigerator functions on DC current. d. Verify the heat exchangers are clean and free of dust and dirt. e. Verify the electrical safety.	The refrigerator cannot function on an AC power supply. The refrigerator cannot function on a DC power supply. Refrigerator does not pass electrical safety tests.